

Rapid, simultaneous analysis of plasma catecholamines and metanephrines by mixed-mode SPE and HILIC LC/MS/MS

Sherri Naughton, Jonathan Danaceau, Erin Chambers, and Kenneth Fountain

Waters Corporation, 5 Technology Dr, Milford, MA 01757

Background: In clinical research, elevated concentrations of plasma catecholamines can be used in conjunction with their O-methylated metabolites (metanephrines) to indicate the presence of conditions such as pheochromocytomas, neuroblastomas, ganglioblastomas and ganglioneuromas. However, these compounds (in particular, norepinephrine, epinephrine, and dopamine) can be a challenge to analyze via reversed-phase LC/MS/MS due to their high polarity. As a result, many research laboratories still analyze this panel using ion-pairing reagents and ECD detection. While reversed-phase LC/MS/MS has been used successfully, challenges still exist due to ion-suppression from matrix components, insufficient retention, and inadequate separation of normetanephrine and epinephrine. This work describes a single extraction and analysis method for monoamine neurotransmitters and metanephrines from human plasma. *Methods:* 250 μ L plasma samples were pretreated with 50 mM $\text{NH}_4\text{CH}_3\text{COO}$, and loaded onto pretreated wells of mixed-mode μ Elution SPE plates. SPE wells were then washed with 20 mM $\text{NH}_4\text{CH}_3\text{COOH}$ and 50:50 ACN:IPA and eluted with 2 x 25 μ L aliquots of 85:15 ACN:H₂O with 2% formic acid. HILIC-based chromatographic separation was achieved using an UHPLC silica-hybrid amide column. MPA and MPB consisted of 30 mM NH_4COO dissolved in 95:5 H₂O: ACN and 15:85 H₂O: ACN, respectively. Compounds were detected by MRM in ESI positive ionization mode. *Results:* All compounds eluted within 2.1 minutes, with baseline separation between normetanephrine and epinephrine enabling their unambiguous identification and quantification. Recoveries ranged from 45-90% and averaged 76%. Matrix effects were less than 25% for dopamine and norepinephrine and under 10% for the remaining analytes. Calibration curves were linear from 10-2000 pg/mL for dopamine, 3-MT, metanephrine and normetanephrine, and from 50-10,000 pg/mL for epinephrine and norepinephrine. Calibration curves for all compounds had R^2 values of 0.999 or greater. %CV and bias values for quality control samples were less than 10% for all analytes at even the lowest QC concentration (40 pg/mL). *Conclusion:* This combination of mixed-mode sample preparation and HILIC chromatography results in a

rapid, robust method with excellent linearity, accuracy, and precision that is suitable for measuring even the lowest endogenous concentrations of these compounds.

For Research Use Only, not for use in diagnostic procedures.